User manual
Benutzerhandbuch
Manual de usuario
Gebruikershandleiding

Microscopes

Article code | Artikelnummer | Código del artículo | Artikelnummer:

760955

BMS S-05-L

WWW.BMSMICROSCOPES.COM


BMS_S-05_GB-DE-ES-NL_manual_v1. PDF9001S05

GB

Thank you for choosing BMS Microscopes.

This stereo microscope enables the study of objects on a larger scale. This basic model operates using a LED battery-operated light source, making it ideal for on-the-go usage.

This highly convenient stereo microscope is suitable for observing insects, minerals or local plants. Developed with the latest advancements in optical and mechanical technology, it is recommended that an adult supervises its use at all times.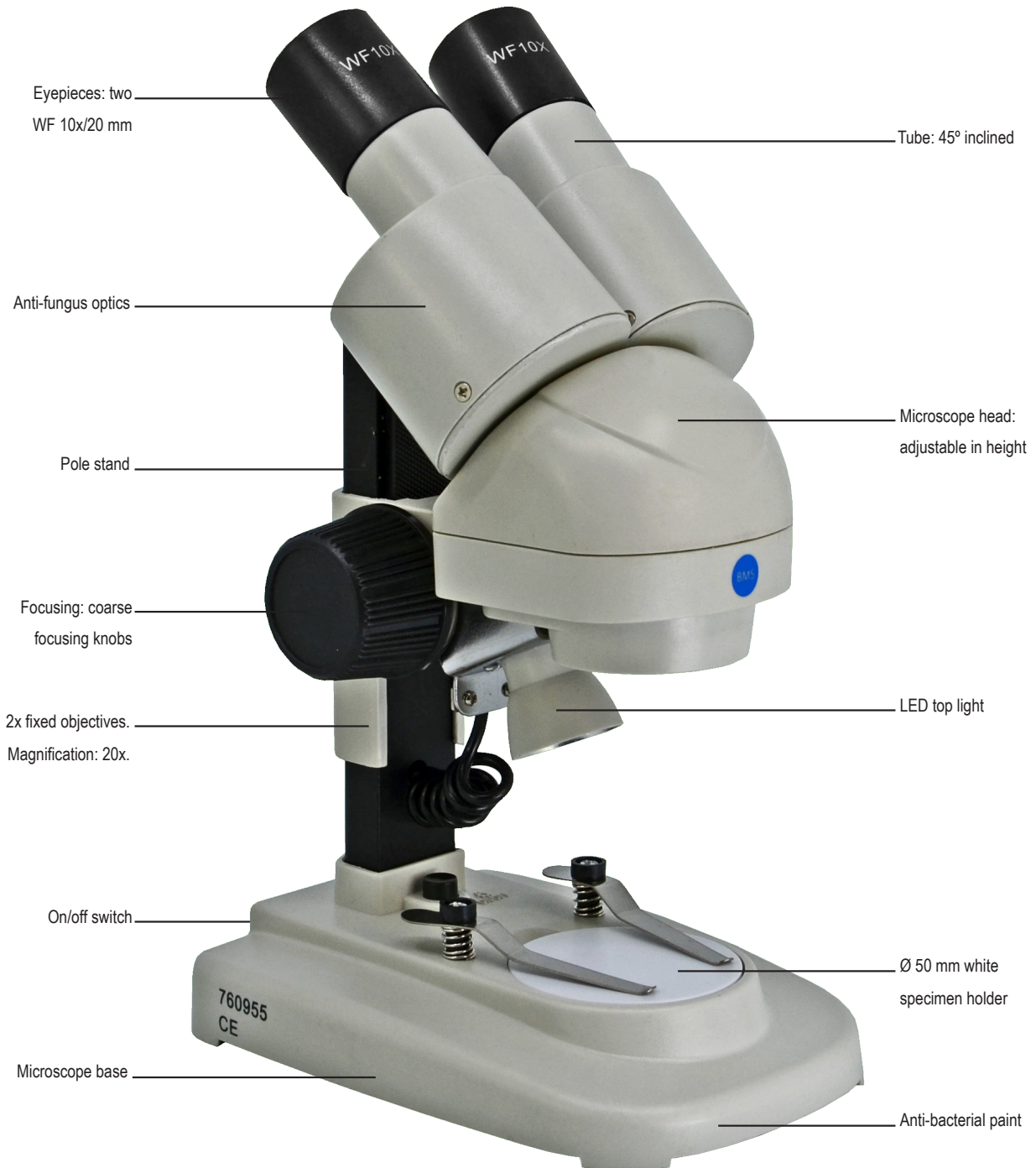

Including: dust cover, manual and Ø 50 mm white specimen holder

Package dimensions (Width x Height x Depth): 17 x 27 x 13 cm

Gross weight: 0,8 kg

Warranty: 5 years

GETTING STARTED

Remove the microscope carefully from the package and retain all packing material. Should you need to return the microscope, you will need the original packing material. Open the battery holder on the bottom of the microscope. Place the batteries in the microscope. Starting with the -side and the +side. Close the battery holder. The microscope is now ready for use.

SAFETY INFORMATION

- Think of the safety of yourself and those being around you.
- Always turn off the microscope before maintenance.
- Work in a clean and proper environment.
- Clean all parts coming in contact with the observed specimen to prevent infection.
- Place the microscope on a sturdy, level table or working surface.
- Always use original accessories.

CAUTION

The delicate components may get damaged if the microscope is not handled with care. Use the microscope in a clean environment. Do not use it when it is subject to direct sunlight, high temperatures, humidity or dust. Carry the microscope with care, do not hold it by the stage. Remove specimen before transport.

CARE & MAINTENANCE

- Microscope is a precision instrument, it should be handled with care.
- The microscope should not be in direct sunlight, but placed in a dry and clean place.
- Avoid high temperature and vibration.
- Avoid dirt and fingerprints on the lens surface, this way your image will not be affected.
- All lenses must be kept clean. Fine dust on the surface of the lens should be blown off with hand blower or wiped off gently with a soft lens tissue. Fingerprints or oil mark on it should be wiped off with a tissue moistened with a small amount of xylene or a mixture of alcohol and ether.
- Never use the organic solution to clean the other surface (especially the plastic surfaces). If necessary, please choose the neutral detergent.
- All lenses are mechanically structured and adjusted to the best state for the microscope, please do not disassemble.
- The rack and pinion should be regularly greased with a small amount.
- After use, cover the microscope with the provided dust cover and store it in a dry and clean place, that is free from moisture to prevent rust.

SPECIFICATIONS


For more specifications, please view the product page on the website:
www.bmsmicroscopes.com

It is within our company policy to develop continuously. BMS Microscopes b.v. reserves the right to change design and specifications without further notice.

BMS Microscopes b.v.
Essebaan 50
2908 LK Capelle aan den IJssel
The Netherlands
Tel.: +31(0)10 458 42 22
E-mail: sales@bmsmicroscopes.com

Environment protection first!

Your appliance contains valuable materials which can be recovered or recycled. Leave it at a local civic waste collection point.

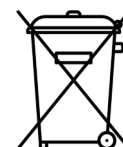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