User manual  
Benutzerhandbuch  
Manual de usuario  
Gebruikershandleiding

Microscopes

Article codes | Artikelnummern | Códigos artículos | Artikelnummers:  
761105, 6762105, 6762535, 6762575, 76250 & 76251

[WWW.BMSMICROSCOPES.COM](http://WWW.BMSMICROSCOPES.COM)


BMS\_stereo-microscopes\_series\_GB-DE-ES-NL\_manual\_v1.PDF9001Stereo

Stereo BMS S-10-2L  
Stereo BMS S-20-2L  
Stereo BMS S-40-2L  
Stereo BMS SL-40  
Stereo BMS ST-50

# GB

Thank you for choosing BMS Microscopes.

Stereo microscopes are widely used in the electronics industry for assembling and testing precision instruments. They are utilized in anatomy, observing and researching animals and plants. They also find applications in schools, scientific research, plant morphology, carving, geology, archaeology and for family research into object shapes.

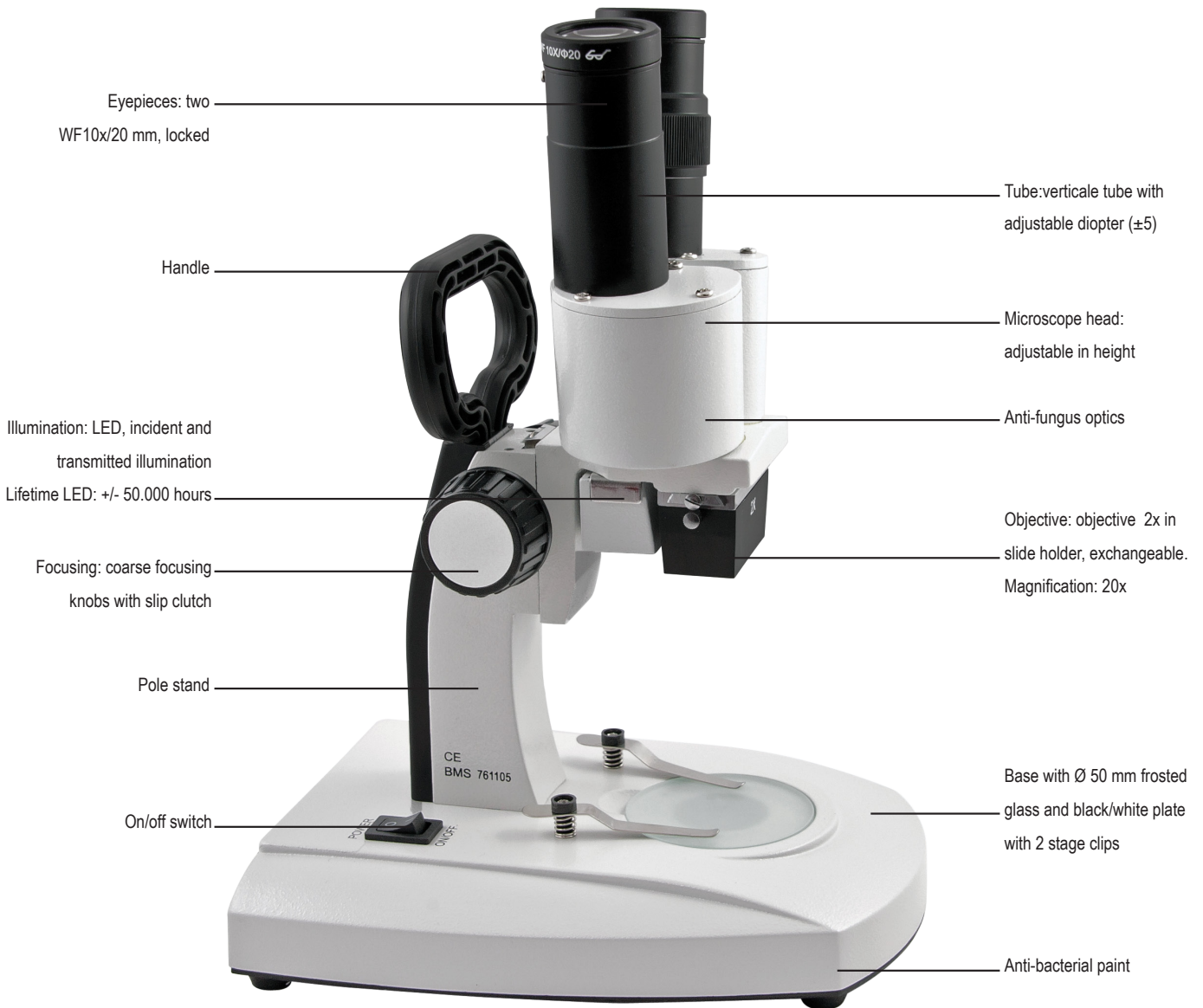Including: dust cover, eyecaps, tension wrench and user manual

Package dimensions (Width x Height x Depth): 29 x 34 x 21 cm

Gross weight: 2,2 kg

Warranty: 5 years

Article code: 761105

Stereo BMS S-20-2L LED (2x) Fixed  
Article code: 6762105


Stereo BMS S-40-2L LED (1x/3x)  
Article code: 6762535

Stereo BMS S-40-2L LED (2x/4x)  
Article code: 6762575

Including: dust cover, eyecaps, tension wrench and user manual  
 Package dimensions (Width x Height x Depth): 29 x 34 x 21 cm  
 Gross weight: 2,6 kg  
 Warranty: 5 years

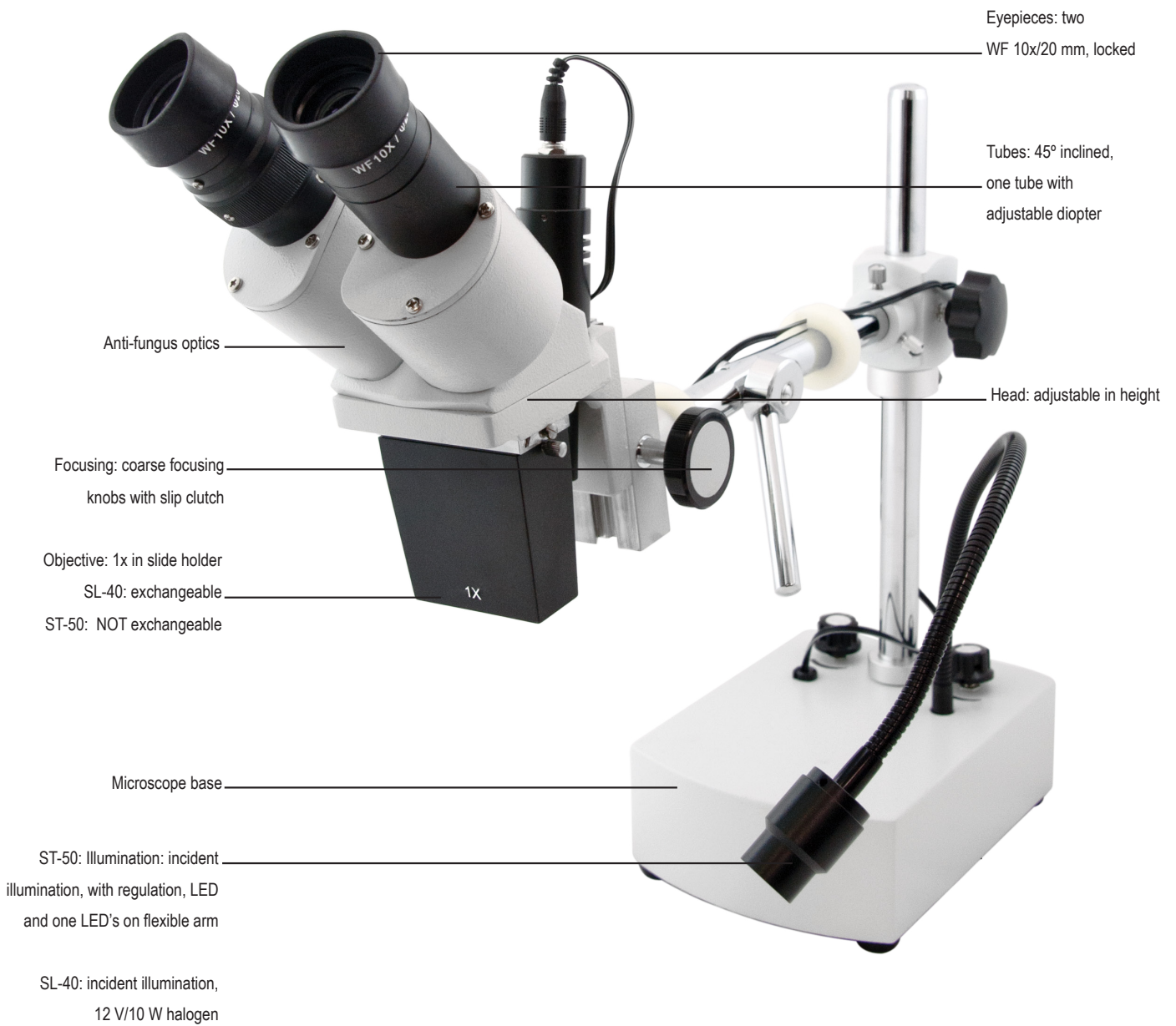Including: dust cover, eyecaps, tension wrench and user manual

Package dimensions (Width x Height x Depth): 41 x 44 x 20 cm

Gross weigh: 6,6 kg (76250 ) and 7,1 kg (76251)

Warranty: 5 years

Article codes: 76250 and 76251

## GETTING STARTED

- Remove the microscope from the packing material.
- Remove and retain all packing material. Should you need to return the microscope, you will need the original packing material.
- Unwrap eyepieces. If included, mount the eye-caps on to the outside of eyepieces.

## MAGNIFICATION

Our Stereo Microscopes offer different magnification options to meet your needs. Depending on the specific type you've purchased:

- Clear 10x magnification (article codes 76250 and 76251).
- Detailed 20x magnification (article codes 761105 and 6762105).
- Versatile 10x/30x (1x/3x) options (article code 6762535)
- Enhanced 20x/40x (2x/4x) magnification (article code 6762575)

## LIGHT CONTROL

- Put 100...240V~ cable into the power socket, open the switch for top/bottom illumination, use the buttons of the illumination and transmission light.
- The frosted glass plate can directly be mounted onto the base, press down gently until secure. To observe the transparent specimen. Use the transmitted light.
- The black and white plate is supplied as an additional item. When using the black and white plate, firstly remove the frosted glass plate and then place the black and white plate over the illumination and press down gently until secure.
- Always put the white side upwards, but if the specimen is white or another bright color, please put the black side upwards as a background to improve the image contrast.

## WHERE TO PLACE THE SPECIMEN

Put the clear specimen in the middle of the working stage, when necessary lock the specimen with the stage clips.

## THE WAY TO USE THE EYE CAP

A pair of eye caps are supplied as an additional item. When using the eye caps, place each eye cap on one eyepiece.

## FOCUS, DIOPTRER ADJUSTMENT, INTERPUPILLARY ADJUSTMENT

The specimen must be placed in the center of the table to be at the correct working position. Then turn the focusing handle slowly until you observe a clear image with the right eyepiece. Then observe the left eyepiece, if the image is not clear, please adjust the ring of the diopter until the image is clear. Adjust both left and right tubes in a way that the distance is appropriate for the eye distance of operator. Once this has been adjusted, you can see a clear image with strong stereo sense.

## POWER OPTIONS

1. Main power cable (100~230V/50Hz), available for EU (item no. 76995), UK (item no. 76999), CH (item no. 76998) and USA (item no. 76996).
2. Rechargeable batteries (AA, 1,5V), only applicable for rechargeable microscopes. It can be ordered as a pack of three (item no. 29554). When this battery pack is placed, it will be charged automatically when the microscope is connected to the mains and the power switch of the microscope is switched on. When the LED indicator is green, the battery pack is fully charged. The microscope can now be used, disconnected from the mains, for 10 hours. When the LED indicator is red, the batteries need to be charged again.
3. Non rechargeable batteries (AA 1,5V, 3 pieces). The microscope will also work with these batteries. **WARNING:** do not connect the microscope to the main electricity when these batteries are placed. This will cause severe damage to the microscope.

## SAFETY INFORMATION

- Always disconnect mains plug from power point before maintenance.
- RISK OF FIRE. Do not charge rechargeable LED microscopes for more than 2 hours.
- Do not disassemble any parts of the microscope, except when noted in the instruction manual. Disassembling the microscope will void the warranty.

- Work in a clean and proper environment.
- Think of the safety of yourself and those working around you.
- Clean all parts coming in contact with the observed specimen to prevent infection.
- Place the microscope on a sturdy, level table or working surface.
- Always use original accessories.

## CAUTION

The delicate components may get damaged if the microscope is not handled with care. Use the microscope in a clean environment. Do not use it when it is subject to direct sunlight, high temperatures, humidity or dust. Carry the microscope with care, do not hold it by the stage. Remove specimen before transport.

## CARE & MAINTENANCE

- Please use our special microscope maintenance kit (item no. 76026).
- Always keep the optical parts clean. Please use the dust cover to protect it.
- Avoid contamination of the objectives.
- In case there is dirt on the lens surface, use distilled water to moist and soft cloth to clean the lens surface. Be very careful and make sure to avoid scratches on a lens surface. Cleaning can also be done by using an alcohol solution or an optical cleaning solution. Ventilate the room. Do not soak the lens in (cleaning) liquid.
- To clean the eyepiece- and objective-lens surface, move the cloth in a spiral motion from the center to the rim.
- Never make a zig-zag movement since this will only spread the dirt.
- To clean painted parts, use also a slightly moistened soft cloth.
- Never use an organic solvent. Use a mild detergent, like soft hand soap.
- Do not remove the objectives from the nosepiece. The objectives were placed in a particular sequence and are adjusted in order to have a parcentered optical system.
- Never dismantle the microscope. This can only be done by factory and/or dealer authorized staff.

## SPECIFICATIONS

For more specifications, please view the product page on the website: [www.bmsmicroscopes.com](http://www.bmsmicroscopes.com)

It is within our company policy to develop continuously. BMS Microscopes b.v. reserves the right to change design and specifications without further notice.

BMS Microscopes b.v.

Essebaan 50

2908 LK Capelle aan den IJssel


The Netherlands

Tel.: +31(0)10 458 42 22

E-mail: [sales@bmsmicroscopes.com](mailto:sales@bmsmicroscopes.com)

*Environment protection first!*

*Your appliance contains valuable materials which can be recovered or recycled. Leave it at a local civic waste collection point.*

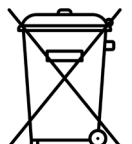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